

—A helping hand for your research

Product Manual

Configuration-specific Monoclonal Antibody Based Cdc42 Activation Assay Kit

Catalog Number: 80701

20 assays

Table of Content

Product Description	3
Assay Principle	3
Kit Components	3
Storage	4
Materials Needed but Not Supplied	4
Reagent Preparation	4
Sample Preparation	5
Assay Procedure	6
Example of Results	8

FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES

NewEast Biosciences 24 Whitewoods Lane Malvern, PA 19355 Fax: 610-945-2008

Phone: 610-945-2007 Web: www.neweastbio.com E-mail: sale@neweastbio.com

Product Description

Small GTPases are a super-family of cellular signaling regulators. Cdc42 belongs to the Rho sub-family of GTPases that regulate cell motility, cell division, and gene transcription. GTP binding increases the activity of Cdc42, and the hydrolysis of GTP to GDP renders it inactive.

Currently the activation of Cdc42 proteins is assayed with the binding of GTP-bound Cdc42 to the p21-binding domain (PBD) of p21-activated protein kinase (PAK). This method is based on the observation that the active, GTP-bound Cdc42 (and Rac) could bind to the PBD of PAK. However, the reproducibility of this method is poor. This is partially due to the relatively quick hydrolysis of GTP to GDP during the assay procedure, and the low binding affinity of PBD to Cdc42-GTP.

NewEast Biosciences Cdc42 Activation Assay Kit is based on the configuration-specific monoclonal antibody that specifically recognizes Cdc42-GTP, but not Cdc42-GDP. Given the high affinity of monoclonal antibodies to their antigens, the activation assay could be performed in a much shorter time. This assay provides the reliable results with consistent reproducibility.

These anti-Cdc42-GTP monoclonal antibody can also be used to monitor the activation of Cdc42 in cells and in tissues by immunohistochemistry.

NewEast Biosciences Cdc42 Activation Assay Kit provides a simple and fast method to monitor the activation of Cdc42. Each kit provides sufficient quantities to perform 20 assays.

Assay Principle

NewEast Biosciences Cdc42 Activation Assay Kit bases on the configuration-specific anti-Cdc42-GTP monoclonal antibody to measure the active Cdc42-GTP levels, either from cell extracts or from in vitro GTPγS loading Cdc42 activation assays. Briefly, anti-active Cdc42 mouse monoclonal antibody will be incubated with cell lysates containing Cdc42-GTP. The bound active Cdc42 will then be pulled down by protein A/G agarose. The precipitated active Cdc42 will be detected by immunoblot analysis using anti-Cdc42 rabbit polyclonal antibody.

Kit Components

1. Anti-active Cdc42, Mouse Monoclonal Antibody (Catalog No. 26905): One vial – 22 μL (1

3

NewEast Biosciences 24 Whitewoods Lane Malvern, PA 19355 Fax: 610-945-2008

Phone: 610-945-2007 Web: www.neweastbio.com E-mail: sale@neweastbio.com mg/ml) in PBS, pH 7.4, containing 50% glycerol and 0.05% sodium azide. This antibody specifically recognizes Cdc42-GTP from all vertebrates.

- 2. Protein A/G Agarose (Catalog No. 30301): One vial 400 μL of 50% slurry.
- 3. <u>5X Assay/Lysis Buffer (Catalog No. 30303)</u>: One bottle 30 mL of 250 mM Tris-HCl, pH 8, 750 mM NaCl, 50 mM MgCl₂, 5 mM EDTA, 5% Triton X-100.
- 4. Anti-Cdc42, Rabbit Polyclonal Antibody (Catalog No. 21010): One vial 22 μL (1 mg/ml) in PBS, pH 7.4, contained 50% glycerol.
- 5. 100 X GTPγS (Catalog No. 30302): One vial -100 μl at 10 mM, use 5 μL of GTPγS for GTP-labeling of 0.5 mL of cell lysate.
- 6. 100 X GDP (Catalog No. 30304): One vial -100 μl at 100 mM, use 5 μL of GDP for GDP-labeling of 0.5 mL of cell lysate.

Storage

Store all kit components at 4°C until their expiration dates.

Materials Needed but Not Supplied

- Stimulated and non-stimulated cell lysates
- 2. Protease inhibitors
- 3. 4 °C tube rocker or shaker
- 4. 0.5 M EDTA, pH8.0
- 5. 1 M MgCl₂
- 2X reducing SDS-PAGE sample buffer 6.
- 7. Electrophoresis and immunoblotting systems
- 8. Immunoblotting wash buffer such as TBST (10 mM Tris-HCl, pH 7.4, 0.15 M NaCl, 0.05% Tween-20)
- Immunoblotting blocking buffer (TBST containing 5% Non-fat Dry Milk or 3% BSA) 9.
- 10. PVDF or nitrocellulose membrane
- 11. Secondary Antibody
- 12. ECL Detection Reagents

NewEast Biosciences Fax: 610-945-2008 4 Phone: 610-945-2007

> Web: www.neweastbio.com E-mail: sale@neweastbio.com

Reagent Preparation

• 1X Assay/Lysis Buffer: Mix the 5X Stock briefly and dilute to 1X in deionized water. Just prior to usage, add protease inhibitors such as 1 mM PMSF, 10 µg/mL leupeptin, and 10 µg/mL aprotinin.

Sample Preparation

Adherent Cells

- 1. Culture cells (one 10-cm plate, $\sim 10^7$ cells) to approximately 80-90% confluence. Stimulate cells with activator or inhibitor as desired.
- 2. Aspirate the culture media and wash twice with ice-cold PBS.
- 3. Completely remove the final PBS wash and add ice-cold 1X Assay/Lysis Buffer to the cells (0.5 1 mL per 10 cm tissue culture plate).
- 4. Place the culture plates on ice for 10-20 minutes.
- 5. Detach the cells from the plates by scraping with a cell scraper.
- 6. Transfer the lysates to appropriate size tubes and place on ice.
- 7. If nuclear lysis occurs, the cell lysates may become very viscous and difficult to pipette. If this occurs, lysates can be passed through a 27½-gauge syringe needle 3-4 times to shear the genomic DNA.
- 8. Clear the lysates by centrifugation for 10 minutes (12,000 x g at 4 °C).
- 9. Collect the supernatant and store samples (~1-2 mg of total proteins) on ice for immediate use, or snap freeze and store at 70 °C for future use.

Suspension Cells

- 1. Culture cells and stimulate with activator or inhibitor as desired.
- 2. Perform a cell count, and then pellet the cells by centrifugation.
- 3. Aspirate the culture media and wash twice with ice-cold PBS.
- 4. Completely remove the final PBS wash and add ice-cold 1X Assay/Lysis Buffer to the cell pellet $(0.5 1 \text{ mL per } 1 \text{ x } 10^7 \text{cells})$.
- 5. Lyse the cells by repeated pipetting.
- 6. Transfer the lysates to appropriate size tubes and place on ice.
- 7. If nuclear lysis occurs, the cell lysates may become very viscous and difficult to pipette. If this occurs, lysates can be passed through a 27½-gauge syringe needle 3-4 times to shear the

5

NewEast Biosciences 24 Whitewoods Lane Malvern, PA 19355 Fax: 610-945-2008

genomic DNA.

- 8. Clear the lysates by centrifugation for 10 minutes (12,000 x g at 4 °C).
- 9. Collect the supernatant and store samples on ice for immediate use, or snap freeze and store at 70 °C for future use.

In vitro GTPyS/GDP Protein Loading for positive and negative controls

Note: In vivo stimulation of cells will activate approximately 10% of the available Cdc42, whereas in vitro GTPγS protein loading will activate nearly 90% of the Cdc42.

- 1, Aliquot 0.5 ml of each cell extract to two microfuge tubes (or use 1 μg of purified Cdc42 protein).
- 2, To each tube, add 20 µl of 0.5 M EDTA (to 20 mM final concentration).
- 3, Add 5 μl of 100 X GTPγS (to 100 μM, final concentration) to one tube (positive control).
- 4, Add 5 µl of 100 X GDP (to 1 mM, final concentration) to the second tube (negative control).
- 5. Incubate the tubes at 30°C for 30 minutes with agitation.
- 6, Stop loading by placing the tubes on ice and adding $32.5 \mu l$ of 1 M MgCl₂ (to 60 mM, final concentration).

Assay Procedure

I. Active Cdc42 Pull-Down Assay

- 1. Aliquot 0.5 1 mL of cell lysate to a microcentrifuge tube.
- 2. Adjust the volume of each sample to 1 mL with 1X Assay/Lysis Buffer.
- 3. Add 1 µl anti-active Cdc42 monoclonal antibody to the tube.
- 4. Thoroughly resuspend the protein A/G Agarose bead slurry by vortexing or titurating.
- 5. Quickly add 20 μL of resuspended bead slurry to each tube.
- 6. Incubate the tubes at 4 °C for 1 hour with gentle agitation.
- 7. Pellet the beads by centrifugation for 1 min at 5,000 x g.
- 8. Aspirate and discard the supernatant, making sure not to disturb/remove the bead pellet.
- 9. Wash the bead 3 times with 0.5 mL of 1X Assay/Lysis Buffer, centrifuging and aspirating each time.
- 10. After the last wash, pellet the beads and carefully remove all the supernatant.

Fax: 610-945-2008 Phone: 610-945-2007

Web: www.neweastbio.com
E-mail: sale@neweastbio.com

- 11. Resuspend the bead pellet in 20 µL of 2X reducing SDS-PAGE sample buffer.
- 12. Boil each sample for 5 minutes.
- 13. Centrifuge each sample for 10 seconds at 5,000 x g.

II. Electrophoresis and Transfer


- 1. Load 15 μ L/well of pull-down supernatant to a polyacrylamide gel (17%). Also, it's recommended to include a pre-stained MW standard (as an indicator of a successful transfer in step 3).
- 2. Perform SDS-PAGE following the manufacturer's instructions.
- 3. Transfer the gel proteins to a PVDF or nitrocellulose membrane following the manufacturer's instructions

III. Immunoblotting and Detection (all steps are at room temperature, with agitation)

- 1. Following the electroblotting step, immerse the PVDF membrane in 100% Methanol for 15 seconds, and then allow it to dry at room temperature for 5 minutes.
 - Note: If Nitrocellulose is used instead of PVDF, this step should be skipped.
- 2. Block the membrane with 5% non-fat dry milk or 3% BSA in TBST for 1 hr at room temperature with constant agitation.
 - Incubate the membrane with anti-Cdc42 polyclonal antibody, freshly diluted 1:50~1000 (depending on the amount of Cdc42 proteins in your samples) in 5% non-fat dry milk or 3% BSA/TBST, for 1-2 hr at room temperature with constant agitation or at 4°C overnight.
- 3. Wash the blotted membrane three times with TBST, 5 minutes each time.
- 4. Incubate the membrane with a secondary antibody (e.g. Goat Anti-Rabbit IgG, HRP-conjugate), freshly diluted 1:1000 in 5% non-fat dry milk or 3% BSA/TBST, for 1 hr at room temperature with constant agitation.
- 5. Wash the blotted membrane three times with TBST, 5 minutes each time.
- 6. Use the detection method of your choice such as ECL.

Example of Results

The following figure demonstrates typical results seen with NewEast Biosciences Cdc42 Activation Assay Kit. One should use the data below for reference only.

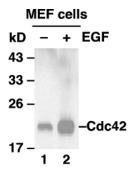

IP: anti-active Cdc42 mAb WB: anti-Cdc42 pAb



WB: anti-Cdc42 pAb

Cdc42 activation assay. MEF cells were treated with (lane 2) or without (lane 1) EGF. Cell lysates were incubated with an anti-active Cdc42 monoclonal antibody (Cat. # 26905) (top panel). The precipitated active Cdc42 was immunoblotted with an anti-Cdc42 rabbit polyclonal antibody (Cat # 21010). The bottom panel shows the Western blot with anti-Cdc42 of the cell lysates used (5% of that used in the top panel).

Phone: 610-945-2007 Web: www.neweastbio.com E-mail: sale@neweastbio.com